


Sanyo MLS-3751L Autoclave

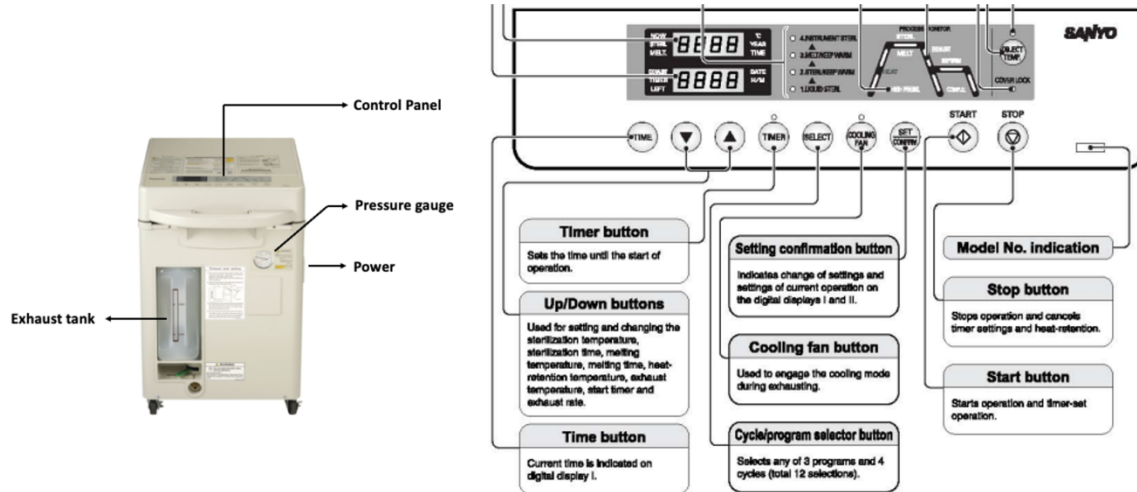

General Operation

1. Fill the exhaust tank with water up to the “Low” level and set the tank to the unit.
2. Turn on the power switch, open the cover.
3. Fill the unit chamber with water until the end of the water level metal of the heater cover is immersed in the water.
4. Put objects to be sterilized in the attached stainless-steel basket, set it in the chamber and closer the cover.
5. After checking the detail of the cycle and settings, press the **START** button.
6. Open the cover and remove the objects after “**Completion**” flickers

Potential Hazards



Preventions to reduce exposure to hazards

